
u-Slide 8 Well high Grid-500

Instructions

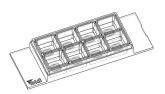

The ibidi product family is comprised of a variety of μ -Slides and μ -Dishes, which have all been designed for high-end microscopic analysis of fixed or living cells. The high optical quality of the material is similar to that of glass, so you can perform all kinds of fluorescence experiments with uncompromised resolution and choice of wavelength. The μ -Slide 8 Well high Grid-500 is an 8 well chamber slide in which cells can be cultivated and, subsequently, investigated with microscopical methods. This open μ -Slide (chambered coverslip) is intended for

cell culture, immunofluorescence, live cell imaging, and high-end microscopy. The Grid-500 is a grid structure for relocating events. It provides 100 distinguishable observation squares of 500 μ m edge length. The grid is clearly visible by phase contrast microscopy and based on the high quality ibidi Polymer Coverslip. The outer dimensions and other parameters are identical to the ibidi μ -Slide 8 Well high.

Material

ibidi μ -Slides, μ -Dishes, and μ -Plates are made of a polymer that has the highest optical quality. The polymer coverslip on the bottom exhibits extremely low birefringence and autofluorescence, similar to that of glass. Also, it is not possible to detach the bottom from the upper part. The μ -Slides, μ -Dishes, and μ -Plates are intended for one-time use and are not autoclavable, since they are only temperature-stable up to 80°C/175°F. Please note that gas exchange between the medium and the incubator's atmosphere occurs partially through the polymer coverslip, which should not be covered.


Optical Properties ibidi Polymer Coverslip

Refractive index n _D (589 nm)	1.52
Abbe number	56
Thickness	No. 1.5 (180 μm)
Material	Polymer coverslip

Please note! The ibidi Polymer Coverslip is compatible with certain types of immersion oil only. A list of suitable oils can be found on page 3.

Geometry

The μ -Slide 8 Well ^{high} Grid-500 provides a standard slide format according to ISO 8037/1.

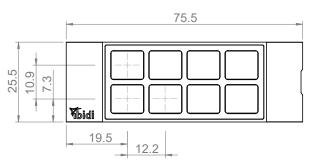

Geometry of the $\mu\text{-Slide8Well}^{high}$ Grid-500		
Outer dimensions in mm $(w \times l)$	25.5 × 75.5	
Number of wells	8	
Dimensions of wells in mm $(w \times l \times h)$	$9.4\times10.7\times9.3$	
Volume per well	300 µl	
Height with/without lid	10.8/9.5 mm	
Growth area per well	1.0 cm^2	
Coating area per well	$2.2\mathrm{cm}^2$	
Bottom	ibidi Polymer Coverslip	

Shipping and Storage

The μ -Slides, μ -Dishes and μ -Plates are sterilized and welded in a gas-permeable packaging. The shelf life under proper storage conditions (in a dry place, no direct sunlight) is listed in the following table.


Conditions	
Shipping conditions Storage conditions	Ambient RT (15–25°C)
	Shelf Life
ibiTreat	36 months

Characteristics of the Grid

The Grid-500 is made of small dots inside the polymer bottom surface of the ibidi $\mu\text{-Slide 8 Well}^{\ high}.$ The structure is imprinted on the side on which cells are growing and does not effect cell growth, coating protocols, or surface properties. Proliferation and cell behavior is comparable with standard non–gridded slides. Cells and grid are in one focal plane.




The grid is made of dots forming grooves, which are $5\,\mu m$ ($\pm\,1\,\mu m$) wide and approximately $5\,\mu m$ deep. Cells can grow in the grooves as well. We recommend using objective lenses up to $20\times$. Anyhow, the optical quality meets the requirements of $63\times$ and $100\times$ oil objective lenses as well (ibidi Polymer Coverslip, No. 1.5).

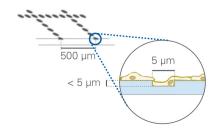

Geometry of the Grid-500

The squares are centred in the wells and indicated by letters and numbers ranging from A to K (J not used) and 1 to 10.

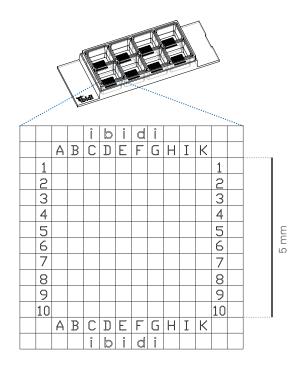

Geometry of the Grid-500		
Number of view fields	100	
Repeat distance	500 μm	
Groove width	5 μm (±1 μm)	
Groove depth	$<$ 5 μ m	
	_ ` ' /	

Surface

The tissue culture-treated ibiTreat surface is a physical surface modification and optimized for adhesion of most cell types. ibiTreat is our most recommended surface modification, because most adherent cells grow well on this hydrophilic version of the ibidi Polymer Coverslip, without the need for any additional coating.

Coating

Detailed information about coatings is provided in Application Note 08: Coating protocols for ibidi labware products.

In short, specific coatings are possible following this protocol:

- 1. Prepare your coating solution according to the manufacturer's specifications or reference.
- 2. Apply 300 µl and leave at room temperature for at least 30 minutes.
- 3. Aspirate the solution and wash with the recommended protein dilution buffer.
- 4. The μ-Slide 8 Well high Grid-500 is ready to be used. Optionally let dry at room temperature. Attention, some coating proteins might degenerate when drying!

Seeding Cells

- Trypsinize and count cells as usual. Dilute the cell suspension to the desired concentration. Depending on your cell type, application of a $5-11 \times 10^4$ cells/ml suspension should result in a confluent layer within 2–3 days.
- Apply 300 µl cell suspension into each well of the slide. Avoid shaking as this will result in inhomogeneous distribution of the cells.
- Cover the slide with the supplied lid. Incubate at 37°C and 5 % CO₂ as usual.

Undemanding cells can be left in their seeding medium for up to three days and grow to confluence there. However, best results might be achieved when the medium is changed every 1–2 days. Carefully aspirate the old medium and replace it by $300\,\mu l$ fresh medium per well.



u-Slide 8 Well high Grid-500

Instructions

Microscopy

To analyze your cells, no special preparations are necessary. Cells can be directly observed live or fixed, preferably on an inverted microscope. The bottom cannot be removed. For optimal results in fluorescence microscopy and storage of fixed and stained samples, ibidi provides a mounting medium (50001) optimized for μ -Dishes, μ -Slides, and μ -Plates.

Tip:

As you may know from 96 well plates, the bent meniscus at the air–liquid interphase in small open wells destroys the phase contrast effect of your microscope image. To avoid this problem, we recommend using our channel Slides such as the $\mu\text{-Slides}$ I Luer and $\mu\text{-Slide}$ VI $^{0.4}$ or a Ph+ Slide.

Chemical Compatibility

The following table provides some basic information on the chemical and solvent compatibility of the μ -Slide 8 Well high Grid-500. For a full list of compatible solvents and more information on chemical compatibility, please visit the FAQ section on ibidi.com.

Chemical / Solvent	Compatibility
Methanol	yes
Ethanol	yes
Formaldehyde	yes
Acetone	yes, without lid
Mineral oil	no
Silicone oil	yes
Immersion oil	See Immersion Oil on page 3.

Immersion Oil

When using oil immersion objectives with the ibidi Polymer Coverslip, use only the immersion oils specified in the table below. The use of any non-recommended oil could damage the ibidi Polymer Coverslip. The resulting leakage may harm objectives and microscope components. All immersion oils that are not listed in the table below should be considered as non-compatible.

Company	Product	Ordering No.	Lot Number	Test Date
ibidi	ibidi Immersion Oil	50101	16-12-27	01/2017
Cargille	Type A	16482	100592	01/2017
Cargille	Type HF	16245	92192	01/2017
Carl Roth	Immersion oil	X899.1	414220338	01/2017
Leica	Immersion Liquid	11513859	n.a.	03/2011
Nikon	Immersion Oil F2 30cc	MXA22192	n.a.	01/2020
Nikon	Silicone Immersion Oil 30cc	MXA22179	20191101	01/2020
Olympus	Silicone Immersion Oil	SIL300CS-30CC	N4190800	01/2017
Zeiss	Immersol 518 F	444960	160706	01/2017
Zeiss	Immersol W 2010	444969	101122	04/2012



Instructions

μ -Slide 8 Well high Grid-500

Ordering Information

$\mu\text{-Slide}$ 8 Well high



Cat. No.	Description
80806	μ-Slide 8 Well ^{high} ibiTreat: #1.5 polymer coverslip, tissue culture treated, sterilized
80806-90	μ-Slide 8 Well high ibiTreat, Bulk Pack: #1.5 polymer coverslip, tissue culture treated, sterilized
80802	μ-Slide 8 Well high Collagen IV: #1.5 polymer coverslip, sterilized
80804	μ-Slide 8 Well high Poly-L-Lysine: #1.5 polymer coverslip, sterilized
80801	μ-Slide 8 Well ^{high} Uncoated : #1.5 polymer coverslip, hydrophobic, sterilized
80807	$\mu\text{-Slide 8 Well}^{\ high}$ Glass Bottom: #1.5H (170 μm ± 5 $\mu m)$ D 263 M Schott glass, sterilized
80807-90	μ -Slide 8 Well ^{high} Glass Bottom, Bulk Pack: #1.5H (170 μm ± 5 μm) D 263 M Schott glass, sterilized

μ-Slide 8 Well high Grid-500



Cat. No.	Description
80806-G500	μ -Slide 8 Well high ibiTreat Grid-500: #1.5 polymer coverslip, tissue culture treated, grid repeat distance 500 μm, sterilized

For research use only!

Further information can be found at www.ibidi.com. For questions and suggestions please contact us by e-mail *info@ibidi.de* or by telephone +49 (0)89/520 4617 0.

© ibidi GmbH, Lochhamer Schlag 11, 82166 Gräfelfing, Germany.